


Download Instructions for DataTrace® RF 1.1.8 software

Please follow the steps below to download DataTrace RF 1.1.8 software from Mesa Laboratories, Inc.

1. Click on this link to download DataTrace RF 1.1.8:
http://www.mesalabs.info/datatrace/DTRF_118.exe
2. Select the option to save the “DTRF_118.exe” file to your local hard drive
3. After the “DTRF_118.exe” file has been successfully downloaded, open up My Computer and then browse to the folder on your local hard drive where this file has been stored.
4. “DTRF_118.exe” is a self-extracting zip file. Double-click on it to extract the files.
5. Once the extraction is complete, a new folder named “DTRF” will be created.
6. Double-click on the DTRF folder to browse its contents. In this folder, the following files and folders can be found:

Docs	File Folder	2/23/2010 1:48 PM
DTRF	File Folder	2/23/2010 1:48 PM
DTW	File Folder	2/23/2010 1:48 PM
Sample	File Folder	2/23/2010 1:48 PM
USBDriver	File Folder	2/23/2010 1:48 PM
AdbRdr90_en_US.exe	34,302 KB Application	7/21/2008 4:44 PM
autorun.inf	1 KB Setup Information	3/12/2008 10:49 AM
Launch.exe	916 KB Application	1/5/2010 10:24 AM
MDAC_TYP.EXE	7,935 KB Application	6/26/1998 12:00 AM

7. Double-click the “Launch.exe” file. The DataTrace 1.1.8 Setup welcome window below will appear:

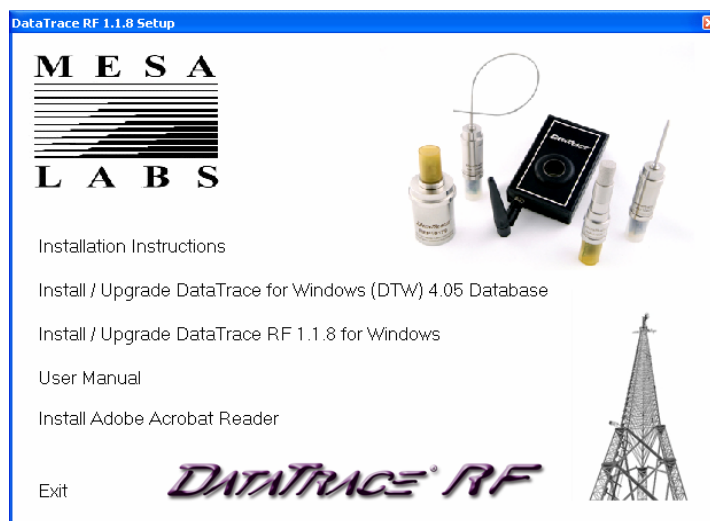

- To install DTRF 1.1.8 on a new computer that does not have any version of DataTrace software installed, please go to page 2 of this document and follow the steps in the **Installation Instructions for DataTrace RF 1.1.8 Software** section.
- To upgrade your existing environment to DTRF 1.1.8, please go to page 3 of this document and follow the steps in the **Upgrade Instructions for DataTrace RF 1.1.8 Software** section.

Installation Instructions for DataTrace® RF 1.1.8 software

These Instructions are applicable to systems that DO NOT have any version of DataTrace Software installed.

To install DataTrace RF 1.1.8 onto a “fresh” system (a computer that does not have any version of DataTrace software installed), please follow the steps below:

1. From the DataTrace 1.1.8 Setup welcome window (shown above), click the **Install/Upgrade DataTrace for Windows (DTW) 4.05 Database** menu option.
2. Then follow the on-screen instructions to install DTW 4.05

Before installing/upgrading DTRF, click the DTW icon on the desktop and verify it opens properly and that a Report can be viewed. If DTW is not functional, do NOT install DTRF. Contact Mesa Laboratories for assistance in debugging a DTW installation issue.

3. Once DTW 4.05 has been successfully installed and verified that it is working properly, click the **Install/Upgrade DataTrace RF 1.1.8 for Windows** menu option on the DataTrace 1.1.8 Setup window.
4. Follow the on-screen instructions to install DTRF 1.1.8
5. Once DTRF 1.1.8 has been successfully installed, click on the **Installation Instructions** menu option on the DTRF 1.1.8 Setup window, then follow the ***Installation Instructions for DataTrace® RF USB Interfaces*** section to install the USB driver for the DataTrace PC Interface.

Upgrade Instructions for DataTrace® RF 1.1.8 software

These Instructions are applicable to systems that already have some version of DataTrace Software installed.

DataTrace RF Software version 1.1.8 requires the DTW database ver. 4.05 be installed and functional prior to the DTRF installation. Access DTW via the desktop icon and determine the DTW version (shown in the upper left corner). If errors occur while accessing DTW, it must be totally uninstalled following the instructions below regardless of the DTW version.

DTW 4.02 or older exists, or, DTW is not functional, or, a Security mode change is desired, Uninstall Required.

If the DTW database contains data that is desired, use the DTW Utilities screen to archive your data. Access your control panel, Add/Remove programs and uninstall / remove DataTrace for windows. When finished, use Windows Explorer and access C:\Program Files\, if a DataTrace folder is found, delete it and all its contents. Access C:\Documents and Settings\All Users\Application Data; if a DataTrace folder is discovered, delete it and all its contents.

If DTRF was present on the computer that had DTW uninstalled, access the Control Panel - Add Remove programs and uninstall DataTrace RF also.

DTW 4.04 exists and is functional, database upgrade required.

Do not uninstall any software. Place the DTRF installation CD in your computer; when the DataTrace RF screen opens, click the second item "Install / Upgrade DataTrace for Windows (DTW) 4.05 Database".

DTW 4.05 exists and is functional, database upgrade required.

Do not uninstall any software. Place the DTRF installation CD in your computer; when the DataTrace RF screen opens, click the second item "Install / Upgrade DataTrace for Windows (DTW) 4.05 Database".

Before installing/upgrading DTRF, click the DTW icon on the desktop and verify it opens properly and that a Report can be viewed. If DTW is not functional, do NOT install DTRF. Contact Mesa Laboratories or your local distributor for assistance in debugging a DTW installation issue.

DTRF 1.1.8 Installation / Upgrade (from DTRF versions older than 1.1.6)

If an old version (1.1.6 or older) of DTRF exists, you will need to uninstall DTRF before proceeding. Use the Add / Remove programs utility to accomplish this if needed. Then install DTRF 1.1.8.

DTRF 1.1.8 Installation / Upgrade (from DTRF 1.1.7 version)

If DTRF 1.1.7 exists, there is no need to un-install DTRF. Place the DTRF installation CD in your computer; when the DataTrace RF screen opens, click the third item "Install / Upgrade DataTrace RF 1.1.8 for Windows".